


## Customize Menu of *Proteus*<sup>®</sup> Analysis

Dr. Alexander Schindler


*Proteus*<sup>®</sup> analysis offers an extensive menu with headlines such as File, Settings, Extras, or Evaluation, which contain a variety of functionalities and options, respectively. As of *Proteus*<sup>®</sup> version 9.0 and higher, a customization of the menu and sub-menus is possible via Settings/Customize Menu... (see figure 1).

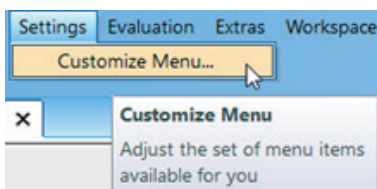

1 Customize Menu in *Proteus*<sup>®</sup> analysis

*Customize Menu* enables to adjust the set of menu items available. In other words, menu items and also the corresponding toolbar icons via checkboxes can be hidden as shown in figure 2. Only menu items checked will be displayed in *Proteus*<sup>®</sup> analysis. Explanations of the settings of this dialogue are available in the help system.

In summary, *Customize Menu* allows for an individualization of *Proteus*<sup>®</sup> analysis according to the users' ideas and requirements.

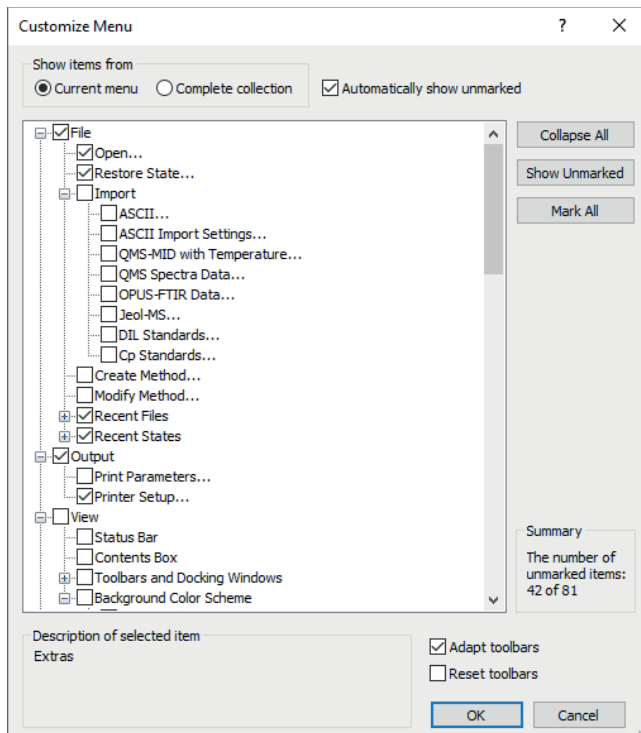

2 Dialog of *Customize Menu* in *Proteus*<sup>®</sup> analysis