APPLICATION NOTE

Investigations of the Influence of Crucible Geometry on the Combustion of Different Carbon Black Samples

Dr. Ekkehard Füglein and Dr. Jan Hanss

Introduction

The methods of thermogravimetry (TGA) are particularly well suited for the investigation of combustion processes. They allow for rapid conclusions regarding the thermal stability of the mostly solid fuel, as well as the reaction temperature and combustion kinetics. Furthermore, both the mass loss during a combustion reaction and the non-combustible mineral ash content can be quantified. Contrary to other reactions such as decomposition or the release of humidity or solvents, combustion is a solid-gas reaction. Therefore, not only must all of the customary parameters such as sample mass, heating rate, and purge gas flow be kept constant, but the measurement results are also influenced by the sample surface, the concentration of oxygen and the crucible geometry, all of which can limit access to the solid sample by the reaction gas.



2 Pierced crucible

Results

For the investigation of carbon black NIST 2975, the crucible types presented in figure 1 were employed. The relationship between the crucible diameter and the filling level of the samples (for the same sample mass) can be seen in figure 3 and table 1.



 $\begin{array}{c} \mbox{I} \quad \mbox{Al}_2 O_3 \mbox{ crucible assortment (from left to right); slip-on plate, short DTA crucible, standard crucible, pierced DTA crucible, mini DTA crucible \\ \end{array}$

To pursue this issue, a series of measurements was carried out with a NETZSCH STA using different crucible geometries under otherwise identical conditions.

The different crucibles are shown in figures 1 and 3; among them is also a pierced DTA crucible which is shown on an enlarged scaling in figure 2 [1].

The investigated carbon black samples are different standards samples such as NIST 2975, Printex 90, activated carbon and carbon balls. These have a diameter of approximately 1 mm to 2 mm and an inorganic structure. The mean particle size of the powder samples is indicated between 20 nm and 50 nm.



3 Crucible assortment, front view (above), top view (below)

Dimensions (mm)	Slip-on plate	Short DTA crucible	DTA crucible	DTA crucible, pierced	Mini DTA*
Ø outer	10	8	8	8	5
Ø inner	10	6	6	6	4

 Tab. 1
 Dimensions of the crucibles shown in in figure 1


*for comparison only; this crucible is not part of the NETZSCH crucible product assortment




APPLICATIONNOTE Investigations of the Influence of Crucible Geometry on the Combustion of Different Carbon Black Samples

When using oxygen as a purge gas, small differences between the various crucible geometries can already be found with regard to the combustion temperature and with regard to the rate of combustion (DTG) (figure 4).


If, however, the concentration of oxygen in the purge gas is reduced to 20% (figure 5) or 5% (figure 6), the crucible geometry appears to play an increasingly important role. The pierced DTA crucible and the slip-on plate obviously allow for a better access of the reaction gas oxygen to the sample. However, the poorer the access of the reaction gas to the solid sample, the greater the tendency for the reaction to shift to higher temperatures and the lower the rate of reaction (DTG). At a nitrogen-tooxygen purge gas ratio of 95:5, the pierced DTA crucible is nearly as "fast" as the slip-on plate. With regard to the reaction behavior, the pierced DTA crucible (figure 2) and the short DTA crucible come closest to the slip-on plate, whereby sample handling for these two crucible types is significantly easier than for the slipon plate.

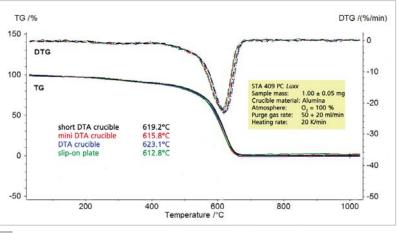



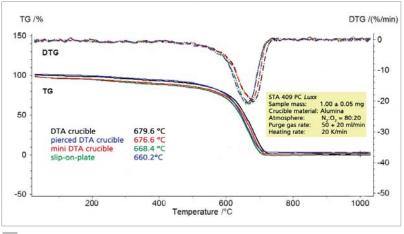

5 TGA-DTG results of NIST 2975 sample (20% O₂)

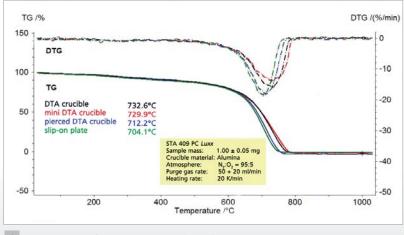


6 TGA-DTG results for NIST 2975 sample (5% O₂)





APPLICATIONNOTE Investigations of the Influence of Crucible Geometry on the Combustion of Different Carbon Black Samples

The dependence of the results on the oxygen content in the purge gas is illustrated in figure 7.


The comparison of different types of carbon black shows significant differences among all of the characteristic values to be determined such as thermal stability, combustion temperature, rate of combustion and residual mass (figures 8 and 9).

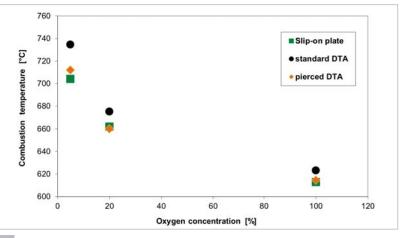

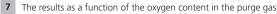


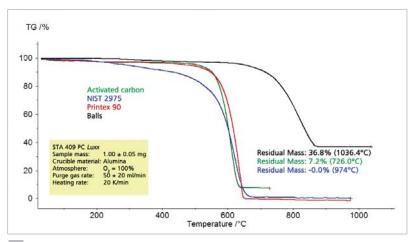

8 Comparison of the four different carbon black samples (TGA, short DTA crucible)

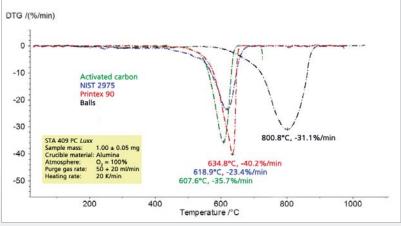

Comparison of the four different carbon black samples (DTG, short DTA crucible)

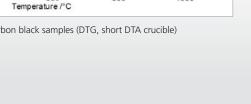

Conclusion

The measurements presented show that the crucible geometry can have significant influence on the interaction between the sample and the purge gas. The combustion reaction of carbon black was used here as an example. Under otherwise identical measurement conditons, as long as the same crucible type was employed within one test series, a comparative evaluation of the samples was possible. The effect of basic measurement conditions, including crucible type, on reaction rate must always be considered when performing kinetic studies. In this case, the slipon plate and pierced crucible proved suitable.

Literature

[1] Protective Anti-Oxidation Coatings for Hot Gas Piping Systems and Their Characterization by Means of a High-Speed Furnace, Thomas Hutsch et.al., NETZSCH OnSet¹⁰, p. 6 – 9.

