

## Coin Cell Cycling in a Novel DSC-Like System

Jean-Francois Mauger, Peter Ralbovsky, Gilles Widawski, and Dr. Peng Ye

### Introduction

Coin cell is a small single cell in the shape of a cylinder or button. It is used in many everyday small portable electronics devices including watches, calculators, remote controls, electronic toys, small LED flashlights etc. It is manufactured in huge quantity every year around the world. There are several different types of coin cells based on chemical composition including silver cell, alkaline cell, zinc-air cell and lithium cell. Some are not rechargeable, while some are rechargeable for example lithium-ion coin cell or Ni-MH coin cell. Compared with the non-rechargeable, disposable cells, the rechargeable cells cost more initially, but the extended lifetime of rechargeable cells make them a better investment for consumers. Plus the rechargeable cells have less impact on the environment as the battery chemicals are often toxic and can pollute environment. Also lithium-ion rechargeable cell offers higher power density and capacity.

Many parameters will influence the performance of rechargeable coin cell including temperature and cycling conditions. It is very important for coin cell manufacturer to know the heat generation during charging/discharging cycles in order to understand the cell energy efficiency and improve cell performance and lifetime. It is also imperative to be able to examine the safety of the whole coin cell when it is subjected to high temperature. This application note shows isothermal cycling of lithium ion coin cell in a novel DSC-like system dedicated to the research and quality control of coin cell battery.

### Instrumentation

The new instrument used is the high temperature coin cell module on NETZSCH MMC 274 *Nexus*<sup>®</sup> (fig 1). It is the only DSC-like system for research and quality control of coin cell battery in the world. The key component of this new instrument is the sensor (fig 2 and 3). It features an innovative differential measurement design based on thermopile for improved sensitivity and stability of heat flow measurement.




1 The high-temperature coin cell module on MMC 274 *Nexus*<sup>®</sup>

It is designed to hold a whole coin cell battery as sample. External battery cyler can be connected to the instrument easily through a plug-and-play LEMO connector. Isothermal cycling of coin cell can be easily performed under various charging/discharging conditions. The new instrument can also do DSC scanning on a whole coin cell which was not possible before.



2 The coin cell sensor assembly

## APPLICATION NOTE Coin Cell Cycling in a Novel DSC-Like System

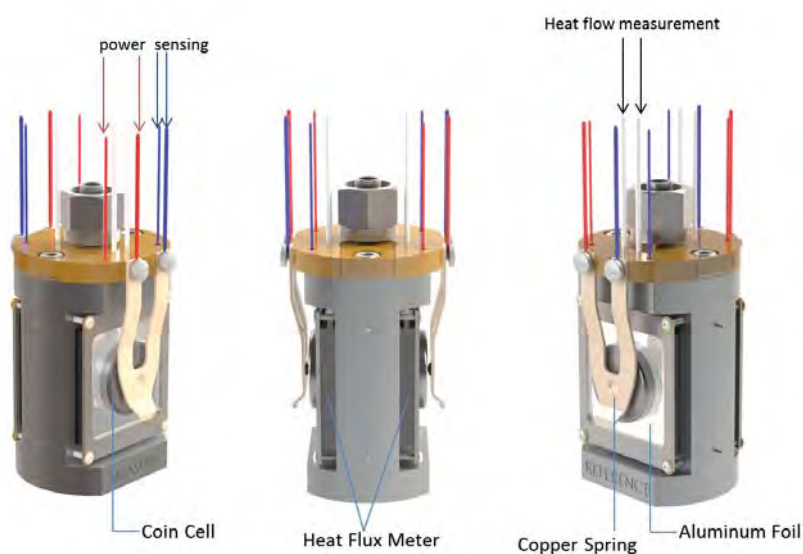


3 Schematic of coin cell sensor

### Results


A LiR2032 lithium ion coin cell battery was used as a sample and an aluminum block of same size was used as reference. The cell was maintained at 40°C throughout the experiment. Constant Current-Constant Voltage (CC-CV) charging and discharging was carried out at 40 mA from 4.2 V to 3.0 V using an external battery cyclers. Several cycles were performed and the heat flow data were quite reproducible. The data shows that charging is character-

ized by a short endotherm followed by a small exothermic reaction while discharging is characterized by a long exothermic reaction (fig. 4). The data can be easily integrated to get the electrical energy in or out of the coin cell and the heat released at the same time (fig. 5). This makes the calculation of battery efficiency straightforward for both charging and discharging processes as shown in figure 4. Without the heat information from HT Coin Cell calorimetry, the efficiency can only be estimated.

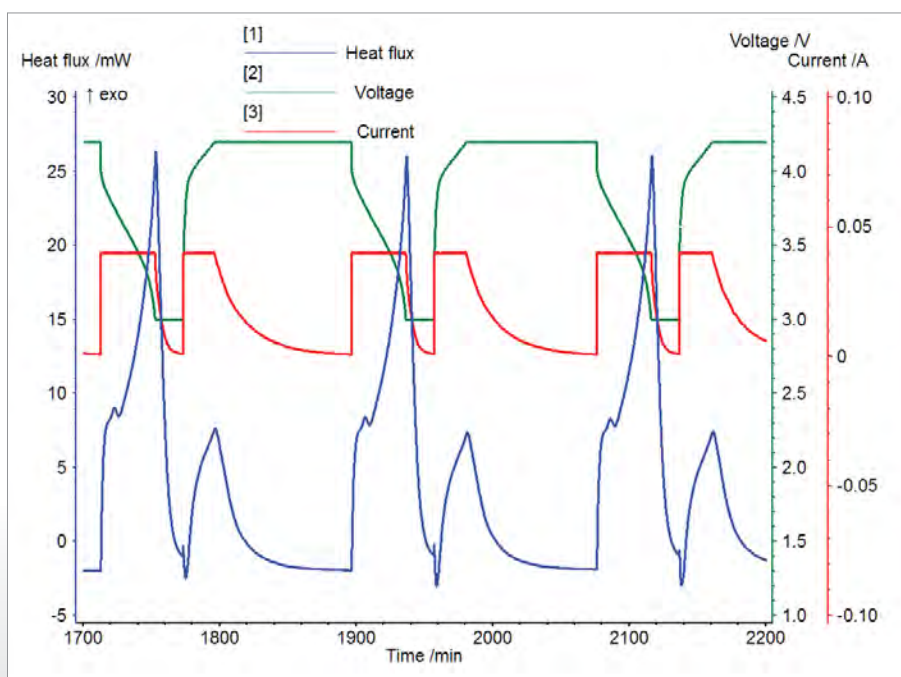

4 Isothermal charging/discharging on coin cell LiR2032 at 40°C

## APPLICATION NOTE Coin Cell Cycling in a Novel DSC-Like System

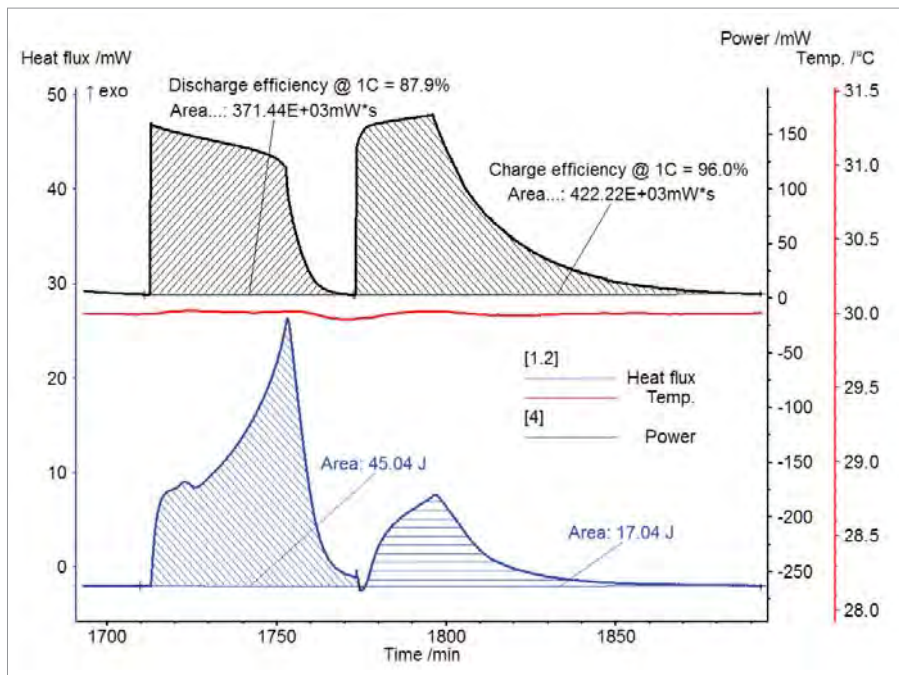

5 Coin cell charging and discharging efficiency calculation using isothermal cycling data

### Conclusion

The new high-temperature coin cell module on MMC 274 Nexus® is a dedicated calorimeter for coin cell analysis. Its innovative sensor incorporates differential principle for high quality data. The sensor is designed to accommodate commercial coin cell battery of different size. A whole coin cell battery can be easily loaded and unloaded. The flexible software enables both isothermal and scanning modes for complete characterization of coin cell.

Thanks to this new instrument, accurate heat flow information can be obtained during charging/discharging test under various conditions. Battery efficiency can be calculated directly with the total heat involved during charging or discharging period. This information is very useful to battery researchers to gain insight into the battery chemistry within the cell and to understand battery performance.